

02019: Programs to run

spdaq20	Program	Command	Comment
	Chip control	<b>gochip</b>	file: 02019_Exp_2.setup
	Pico shaper/discriminator control	<b>gopico</b>	LASSA: input-files/02019/ WU-Shap.dat and WU-Disc.dat
			Target Si: input-files/02019/WU-Target-Shap.dat
			Target Si: input-files/02019/WU-Target-Disc.dat

spdaq29	Program	Command	Comment
	Elog server	<b>goelog</b>	

Any data U6 PC (linux)	Program	Command	Comment
	Readout	<b>godaq</b>	
	Scalers	<b>goscaler</b>	
	Monitoring (V, T...)	<b>goepics</b>	
	Elog client	<b>elog</b>	The server has to be started on spdaq29.
	SpecTcl	<b>gospec</b>	
	Bias for HiRA Si and CsI (CAEN)	<b>telnet 35.9.56.160 1527</b>	
	id = user; passwd = user		
	tab to access menu		
	goup 1 = Si; group 2 = CsI		